



Operation Manual for Spindle Code. L0

1. List for Spindle Code. L0

- 1.1 Joint rod 1
- 1.2 Hook rod 1
- 1.3 Frame 1
- 1.4 Spindle Code. L0 1
- 1.5 Pipe 1

2. Installation

- 2.1 Install the frame on the viscometer;
- 2.2 Connect the joint rod with the hook rod, also the hook rod is in connected with Spindle Code. L0;
- 2.3 Install the joint rod on the joint screw;
- 2.4 Push the measured liquid into the pipe and then fix the frame with the screw , ensure that Spindle Code.0 is covered with the measured liquid;
- 2.5 Set the viscosity meter according to “Operation Manual for NDJ Series and SNB Series Viscometer”, then you can see the viscosity of the measured liquid.

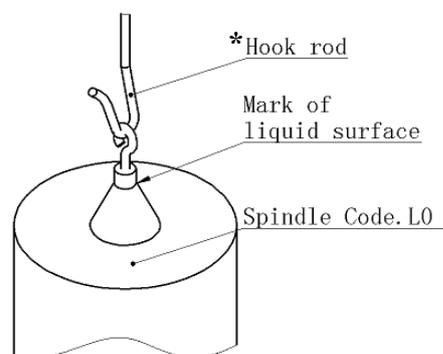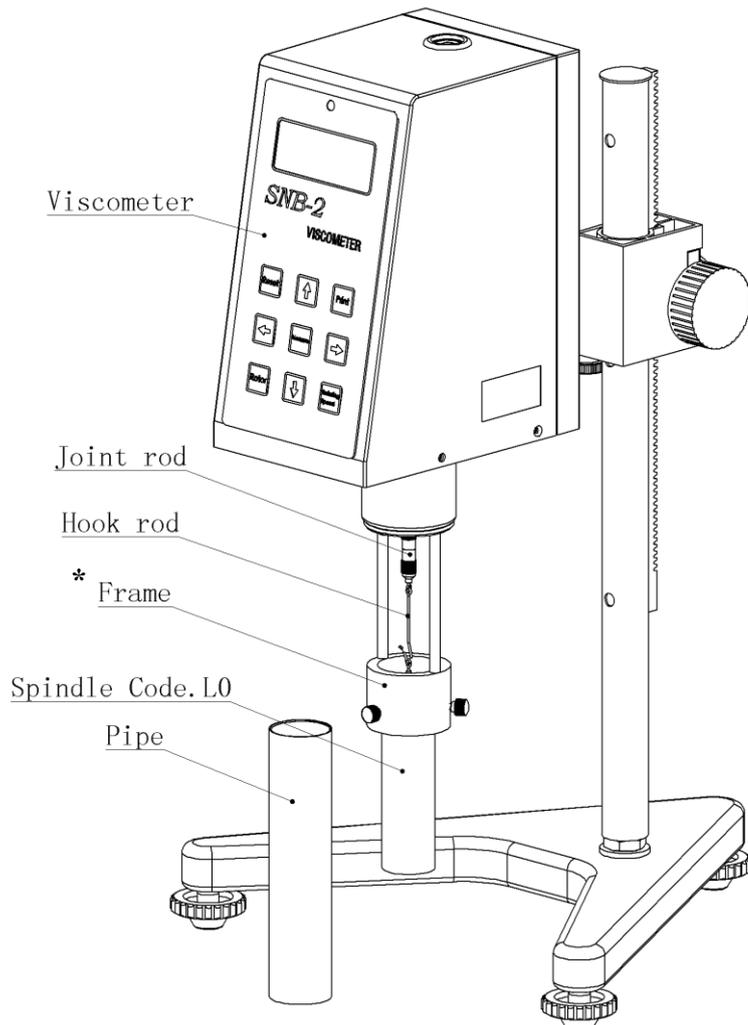

*Note:

- 1. Pay attention to the opening direction of the hook rod and the rotor hook.
- 2. Pay attention to the direction of the screws on the frame (one facing forward).